

Case study - Yeast

Rapid methods to extract DNA & RNA from *Cryptococcus neoformans*.

Alessandro Bolano, Silvia Stinchi, Roberta Preziosi, Francesco Bistoni, Massimo Allegrucci, Franco Baldelli, Alessandro Martini, Gianluigi Cardinali.

FEMS Yeast Research. 2001. Vol 1.

Introduction

Extraction of nucleic acids from the pathogenic yeast *Cryptococcus neoformans* is hampered by a thick and resistant capsule accounting for at least 70% of the whole cellular volume.

This study presents an effective procedure based on mechanical cell breakage using the FastPrep® system to extract RNA from *C. neoformans* and other capsulated species.


Overview

- **Keywords:** Yeast, DNA, RNA, Extraction, Capsule, Method, *Cryptococcus neoformans*
- **Aim of the study:** Development of consistent extraction method for DNA & RNA extraction from resistant strains of *Cryptococcus neoformans*
- **Application:** RNA extraction
- **Sample name:** *Cryptococcus neoformans*
- **Sample type:** Yeast
- **Material:** FastPrep-24™ instrument, 0.4-0.6 mm glass beads
- **Buffers:** TE buffer, RNA lysing solution (0.5% SDS and 0.5% N-laurylsarcosine)

Protocol and Parameters


Cells were grown in YEPD (yeast extract 1%, peptone 1%, dextrose 2%) for 18h at 25 °C in 500-ml shaken (150 rpm) flasks with a liquid to air ration of 1: 10.

1. Cells from 15 ml overnight culture (approx.10⁹) were collected, washed with cold water, resuspended in 200 µl of TE and distributed into two 1.7 ml microcentrifuge tubes.
2. 0.5 ml of glass beads (0.4-0.6 mm), 250 µl of RNA lysing solution, 250 µl 4 M guanidine thiocyanate with 25 mM sodium citrate, pH 7, 0.1 M β-mercaptoethanol, 500 µl phenol, pH 5 and 100µl chloroform/isoamyl alcohol (24 :1) were added to each tube.
3. Both tubes were placed in the FastPrep-24™ instrument and processed in 4 cycles of 40 sec each. Between the cycles samples were placed on ice.
4. After last cycle tubes were removed from the instrument, placed 5 min in ice and spun 10 min at 13,000 x g.
5. The upper phase was transferred to fresh microcentrifuge tube for RNA purification.

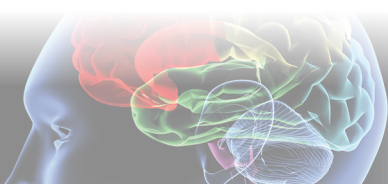

Results

Effective *C. neoformans* RNA extraction from clinical isolates

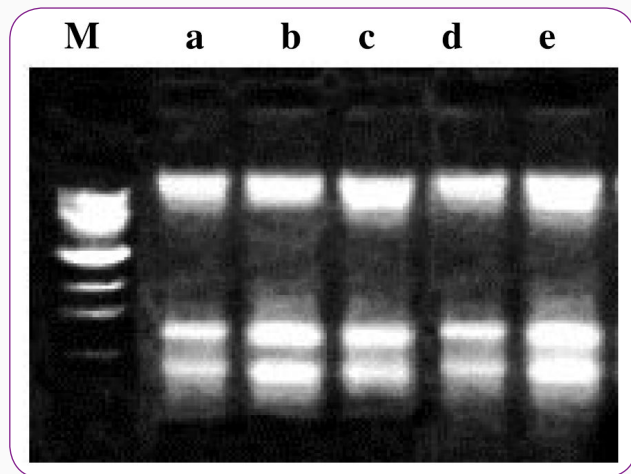


Electrophoresis of total cellular RNA extracted from C. neoformans and S. cerevisiae on a 0.8% agarose.

Lane M, 1 kb ladder. Lanes a, b, c and d, C. neoformans RNA extracted from clinical isolates from cerebrospinal fluid of an AIDS patient (a, first episode; b, relapse) and from CBS collection (c, CBS 6995 encapsulated strain; d, CBS 7698 non-capsulated strain). Lane e, S. cerevisiae RNA extracted from the DBVPG 6820 strain.

Conclusion

- RNA purification is accomplished using FastPrep® system and glass beads after a preliminary bead beating treatment
- Yields range around 1 mg RNA from 15 ml overnight culture (10^9 cells)
- RNA appears undegraded, making it suitable for molecular manipulations

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Successful sample preparation using the MP Biomedicals FastPrep® product line has been highlighted in thousands of scientific articles. To access articles and other materials, visit www.mpbio.com/FastPrepLibrary.



www.mpbio.com



33F0715-1